

LLM-Driven Library Screening for Discovery of Antibody-Based GLP-1R Agonists

Georgia Pentland, Olivier Touzelet, Lorna Stewart, Conor Breen, Sean Tierney, Giulia Mignone, Brendan Clarke, Nicola Hamilton, Simon Larsen, Richard Buick, Adrian Kinkaid.

Fusion Antibodies plc., Belfast, Northern Ireland.



Background

LLM-based Antibody Discovery^[1,2]

- Antibody-based agonists offer potential advantages over peptide agonists, including longer half-life, improved specificity, and favourable biodistribution
- Multi-pass transmembrane GPCRs remain challenging targets for antibody discovery, with relatively few FDA approvals to date. Structurally constrained extracellular epitopes can limit antibody binding and functional modulation of receptor signalling
- Recent advances in large language models (LLMs) have accelerated de novo antibody generation, increasing the potential for successful campaigns

GLP-1R as a Therapeutic Target

- Glucagon-like peptide-1 receptor (GLP-1R), a class B G protein-coupled receptor (GPCR), is a validated therapeutic target for type 2 diabetes and obesity
- Aim of isolating agonistic GLP-1R-binding antibodies capable of activating downstream cAMP signalling to promote glucose-dependent insulin secretion

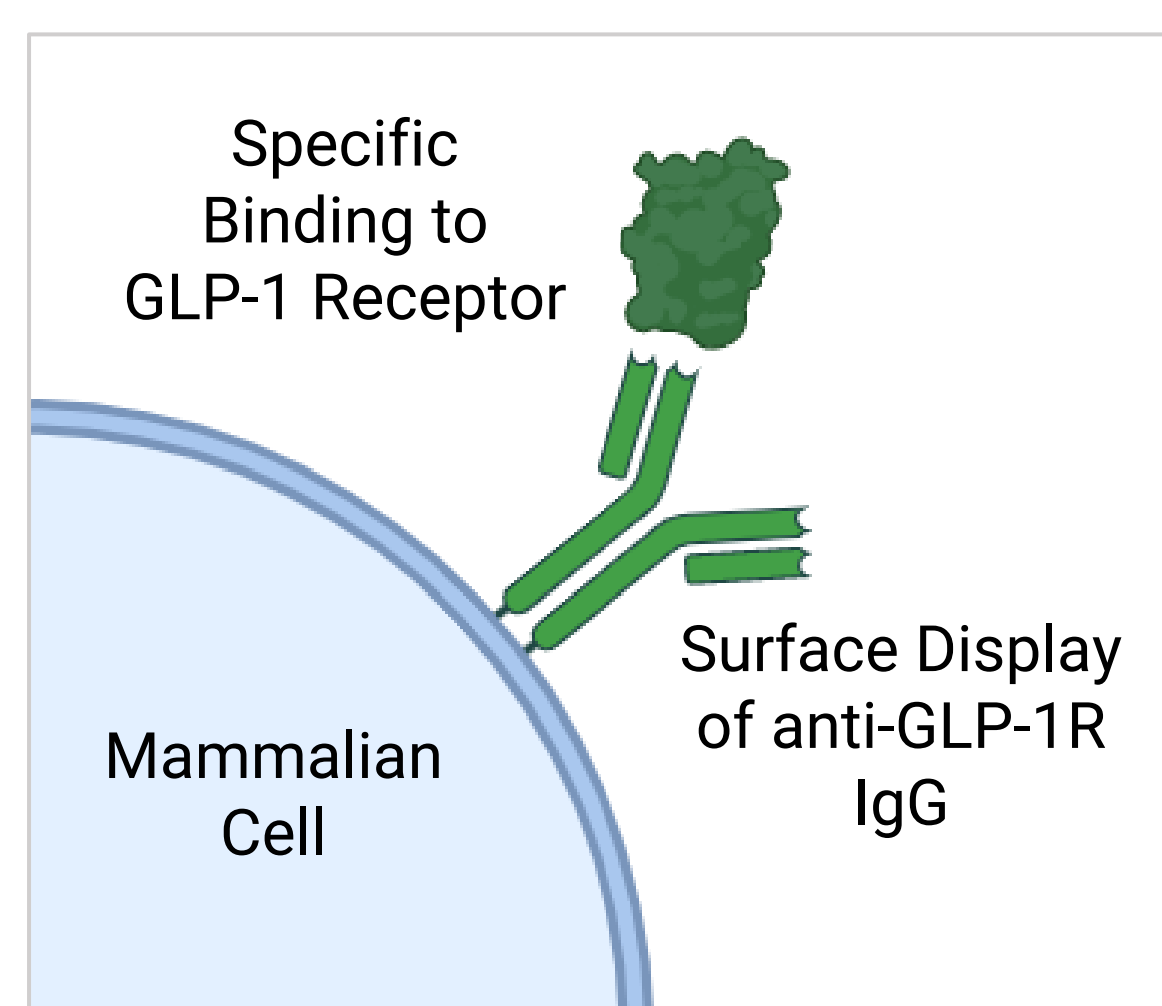


Fig 1. Mammalian cell displaying a full-length anti-GLP-1R IgG that is binding specifically to the GLP-1 Receptor

AI/ML-ab™ In-Silico Design Platform

- Traditional discovery of high affinity antibodies relies on time-consuming immunisation and manual engineering, including CDR grafting and mutation
- AI/ML-ab™ preserves native eukaryotic folding and post-translational modifications whilst facilitating selection for binding, biophysical properties, and reduced immunogenicity
- Combination of AI/ML-ab™ and OptiMAL® platforms enable screening of diverse AI/ML-designed libraries within high-throughput mammalian display

Methodology

LLM Training

- Interrogation of >2 billion human antibody sequences from large-scale databases, including OPIG Observed Antibody Space (OAS), enabling learning from naturally occurring repertoires
- Generation of 3,961 heavy-chain and 1,913 light-chain sequences, combinatorially paired to produce ~8 million unique antibody candidates predicted to bind full-length human GLP-1R
- Sequences were assessed for developability as drug candidates and prioritised according to desired sequence and germline characteristics

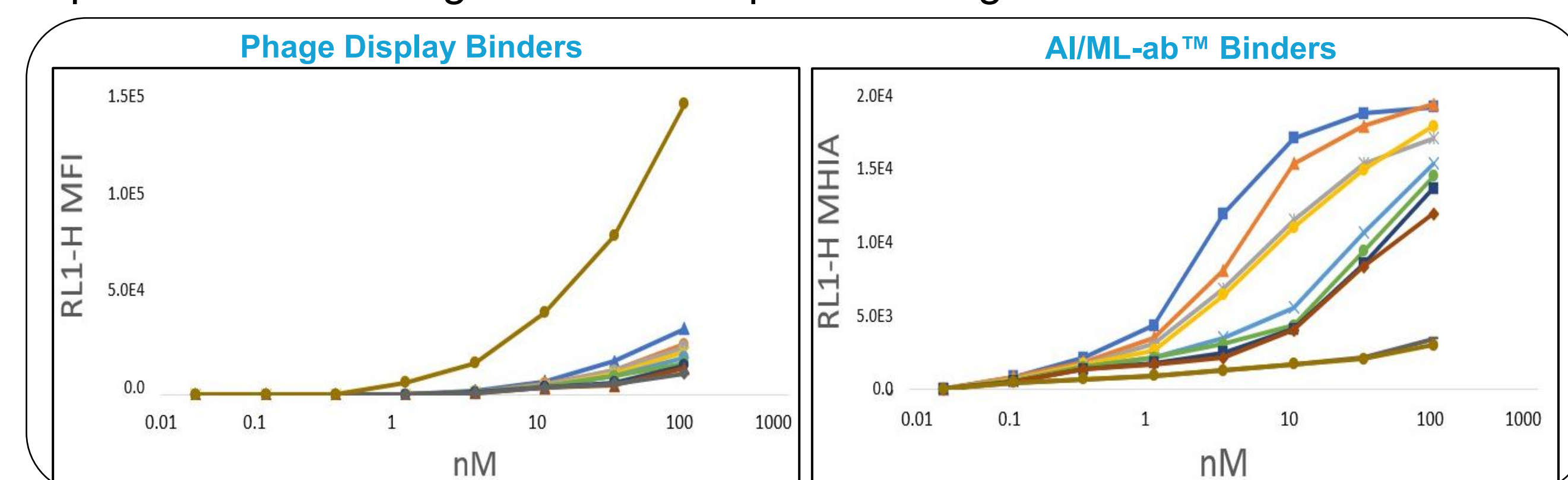


Fig 2. Comparison of Binding Profiles in Phage Display (left) and AI/ML-ab™ in-silico Libraries (right). AI/ML-ab™ binders exhibited greater concentration-dependent affinities.

Library Generation

- Library incorporated into HEK293T via lentiviral transfection and transduction with GFP as a reporter for surface display of anti-GLP-1R IgG variants
- Iterative Magnetic- and Fluorescence- Activated Cell Sorting (MACS) and (FACS) conducted to enrich GLP-1R-binding antibodies, incorporating positive and negative selection with recombinant GLP-1R protein
- Enriched hits were purified and re-expressed as recombinant H + L pairs for biophysical and functional characterisation, assessed against the theoretically-ranked top 96 sequences



Results

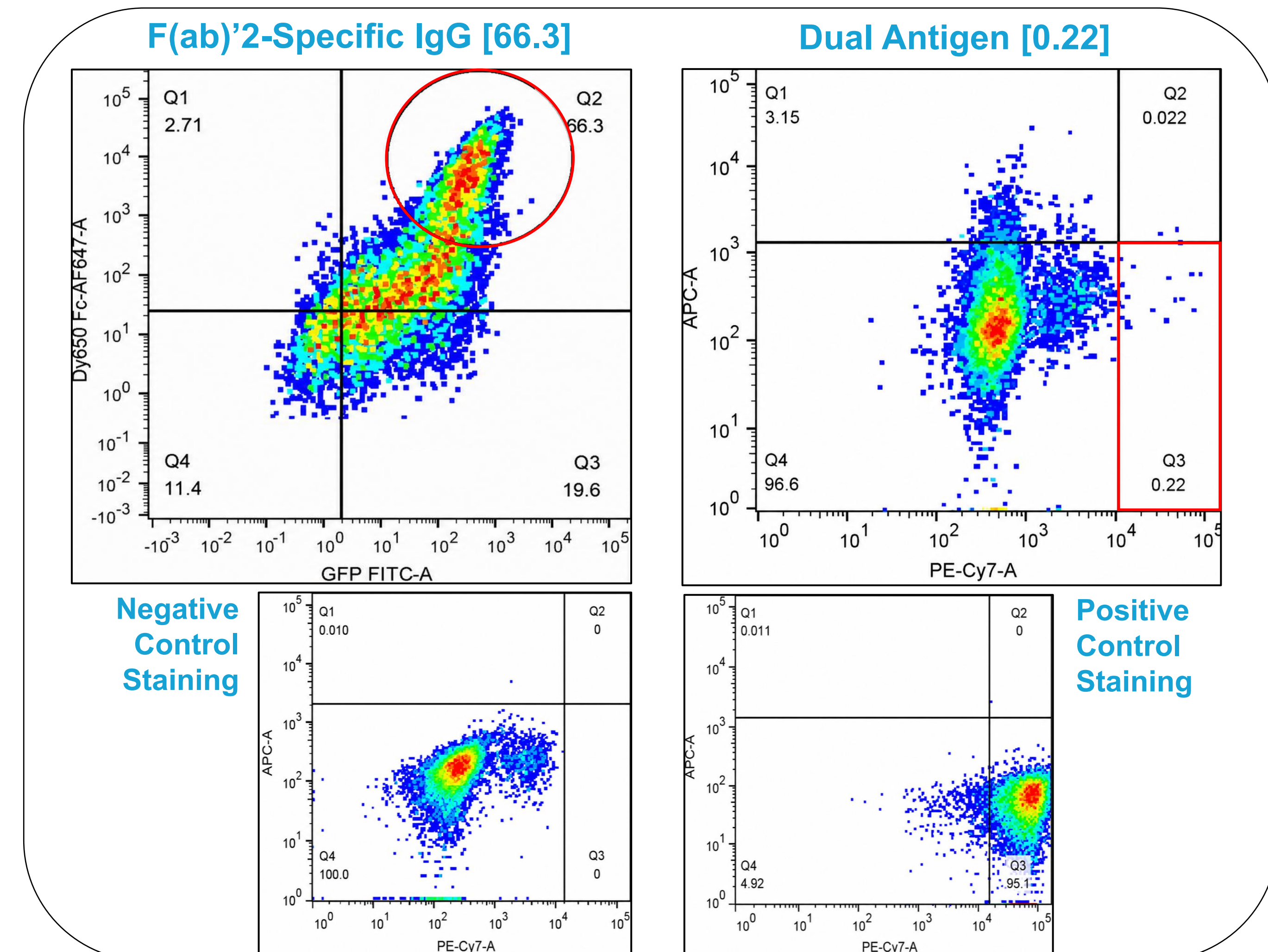


Fig 3. Selection of GLP-1R-specific Binders via FACS. Following iterative antigen-specific enrichment via Magnetic-Activated Cell Sorting (MACS), an enriched population of **GFP positive cells** was detected by flow cytometry. An initial bulk sort to isolate full-length IgG variants (via anti-hu F(ab)'2 IgG Display) was conducted, yielding **F(ab)'2-Specific IgG [Q2, 66.3%]**. Post-recovery, the bulk-sorted population underwent **dual-antigen staining**, using GLP-1R recombinant protein for positive selection (**PE-Cy7-A**) and a negative antigen control (**APC-A**). **GLP-1R-specific binders** were isolated whilst excluding non-specific antigen binding [**Q3, 0.22%**]. Selected events were single-cell sorted (SCS) into 96W plates, expanded, and validated for antigen binding by flow cytometry (FC).

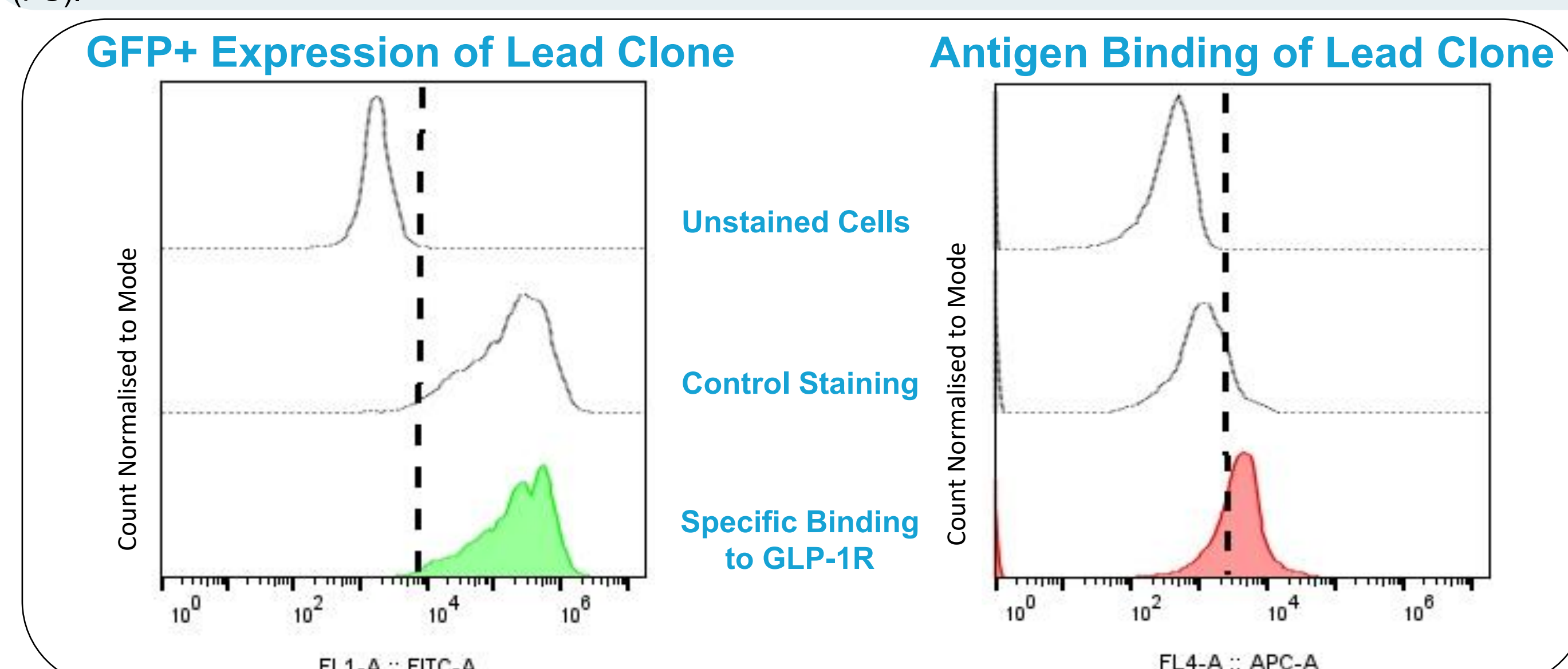


Fig 4. Flow Cytometry Validation of Identified GLP-1R Antibody Hits. Individual clones brought forward from Single Cell Sorting (SCS) were assessed by flow cytometry (FC) for GFP expression and antigen binding to anti-GLP-1R. **GFP+** fluorescence indicated stable presence of the antibody library, whilst **APC+** fluorescence confirmed surface display of antibodies capable of binding to the target recombinant protein (GLP-1R). Identified hits demonstrated **dual GFP+/APC+** signal. Unstained HEK293T cells and a non-targeting antibody stain control showed no detectable APC signal, supporting **GLP-1R specificity**. Rescreened clones were subsequently cryopreserved and re-expressed for downstream characterisation and NGS analysis.

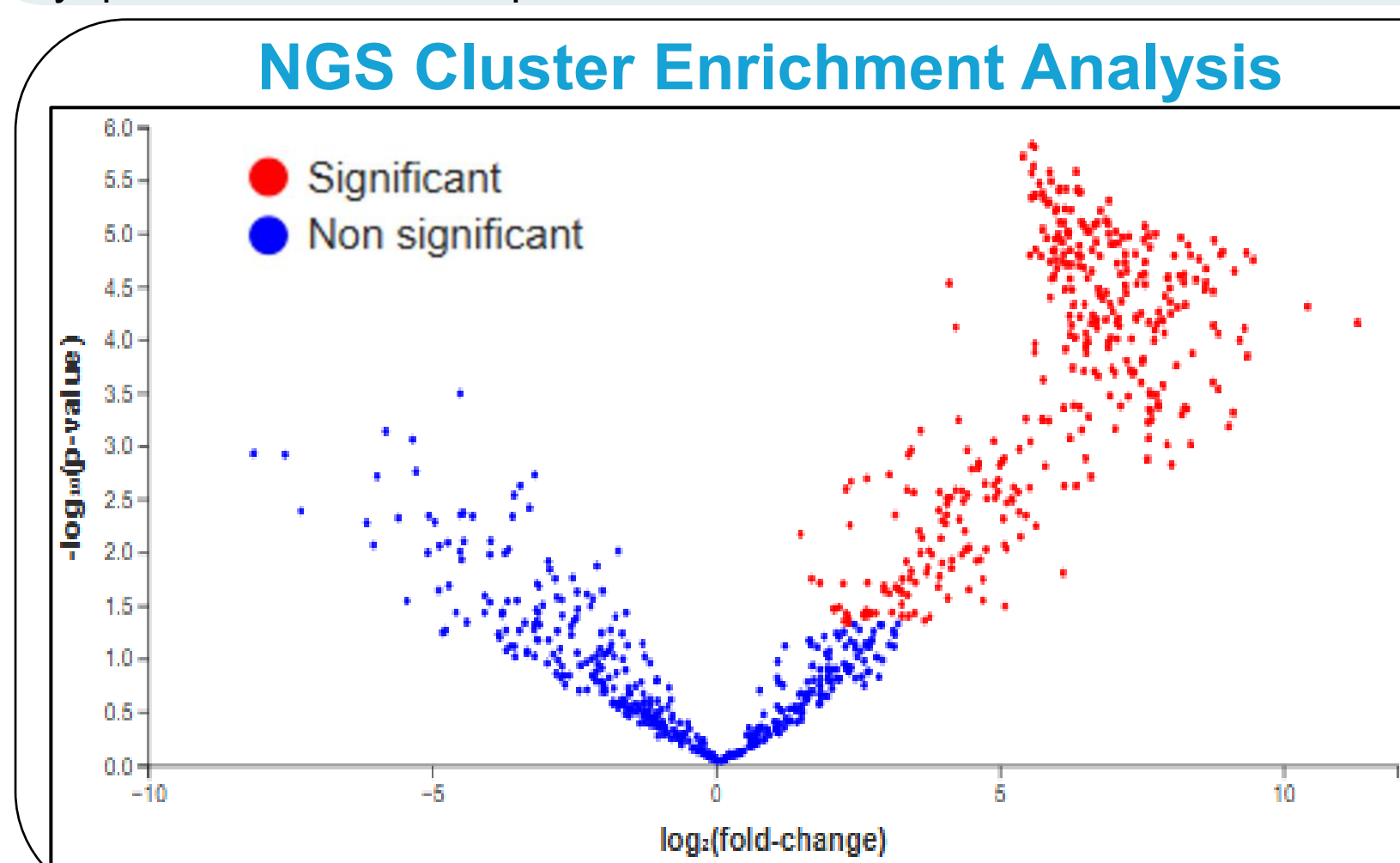


Fig 5. Cluster Analysis of Enriched IgG Sequences. Antibody sequence clusters identified following iterative MACS rounds were compared with the unprocessed library. **Red dots (n=435)** represent **statistically significant positive fold-change**, whilst **blue (n=341)** remain **unenriched or depleted**. Overall enrichment was observed; however, insufficient specificity for the desired GLP-1R binding phenotype remained. Five out of 96 top-ranked hits were contained within the enriched dataset, though only one demonstrated on-target binding.



Conclusions & Future Work

- Combining LLM-driven de novo antibody generation with high-throughput mammalian display is a promising approach to overcome key obstacles associated with discovery against challenging targets
- GLP-1R-binding antibody sequences were successfully recovered from the de novo library, validating the existing methodology and highlighting the potential of AI/ML technologies to shape the future of antibody discovery
- Optimisation of enrichment and selection stringency is necessary to increase target-specific enrichment for GLP-1R. Further characterisation studies will aim to ultimately isolate functional GLP-1R agonists

Acknowledgements Thank you to the team at Fusion Antibodies who have been a part of the development of the AI/ML-ab™ In-Silico Antibody Discovery platform to date. Thank you to Prof. Christopher Scott and the Fusion Antibodies Scientific Advisory Panel for their advice. Figures were created using BioRender, FlowJo, and PipeBio.

References (1) Kim J, McFee M, Fang Q, *et al.* Computational and artificial intelligence-based methods for antibody development. *Trends Pharmacol Sci.* 2023;44(3):175-189. (2) Bai G, Sun C, Guo Z, *et al.* Accelerating antibody discovery and design with artificial intelligence: recent advances and prospects. *Semin Cancer Biol.* 2023;95:13-24.